

ZEISS Axio Zoom.V16 Quick Guide

Initial Set Up

1 Incident Light Source HXP 200 C



2 Transmitted Light (turn the rotary knob clockwise)



3 Electronic Module EMS 3 (turn the device on in the back)



4 Computer




5 Wait for SYCOP 3 to initialize




6 Start ZEN 2.1 (blue edition)



7 Place the specimen into the round insert plate

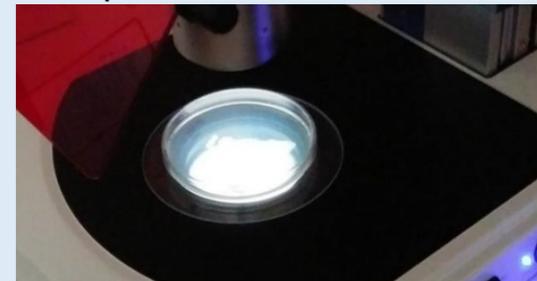


8 Set the tube for camera acquisition

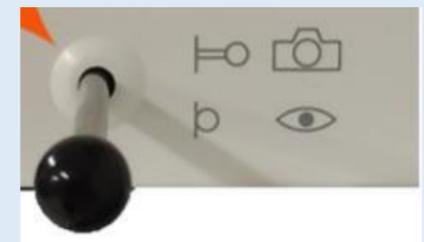

Simple Acquisition


1 In the Locate tab, select the contrast method or filter set

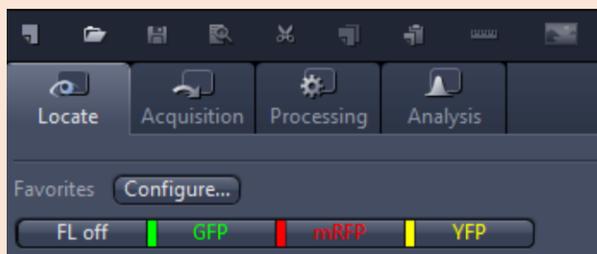

2 Operate the diaphragm slide accordingly



3 Press Live. Set the exposure time

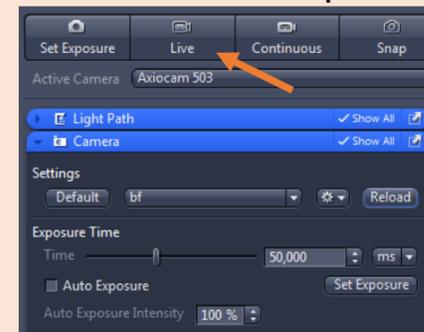


4 Focus the sample using SYCOP 3. Press Snap to acquire



Final Set Up

1 Remove the specimen from the round insert plate

2 Press the Shut Down button on SYCOP 3

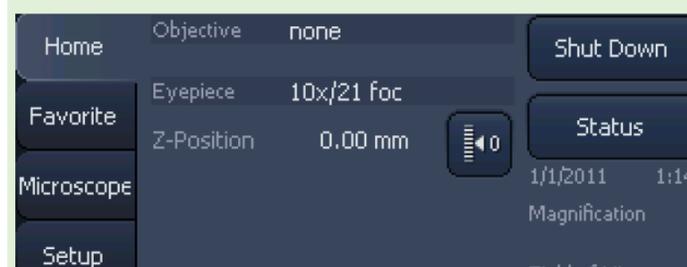

3 Save the data and switch the devices off in the reverse order of the Initial Set Up

4 Sign the logbook before you exit the room